gibco

invitrogen

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Epi5™ Episomal iPSC Reprogramming Vectors

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Product	Size	Cat. No.
CTS™ CytoTune™-iPS 2.1 Sendai Reprogramming Kit	1 Kit	A34546
CytoTune™-iPS 2.0 Sendai Reprogramming Kit	1 pack	A16517
CytoTune™-iPS 2.0 Sendai Reprogramming Kit	3 pack	A16518
Epi5™ Episomal iPSC Reprogramming Kit	1 kit	A15960
CytoTune™ EmGFP Sendai Fluorescence Reporter	1 pack	A16519
Episomal iPSC Reprogramming Vectors	1 tube	A14703

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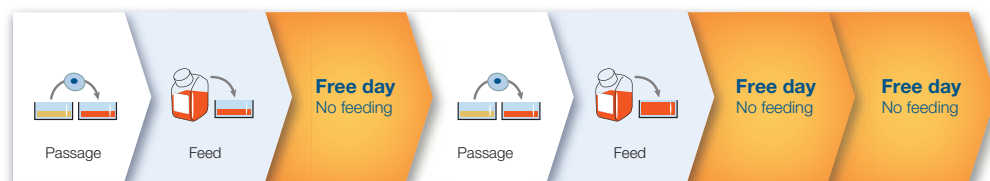
Essential 8™ Medium & Essential 8™ Flex Medium

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The Flex way: Eliminate daily feeding

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Product	Size	Cat. No.
StemPro™ BM Mesenchymal Stem Cells	1 x 10 ⁶ cells	A15652
StemPro™ Human Adipose-Derived Stem Cell Kit	1 kit	R7788110
StemPro™ Human Adipose-Derived Stem Cells	1 vial	R7788115
MesenPRO RS™ Medium	1 kit	12746012
StemPro™ MSC SFM	1 kit	A1033201
StemPro™ MSC SFM XenoFree	1 kit	A1067501
StemPro™ Adipogenesis Differentiation Kit	1 kit	A1007001
StemPro™ Chondrogenesis Differentiation Kit	1 kit	A1007101
StemPro™ Osteogenesis Differentiation Kit	1 kit	A1007201

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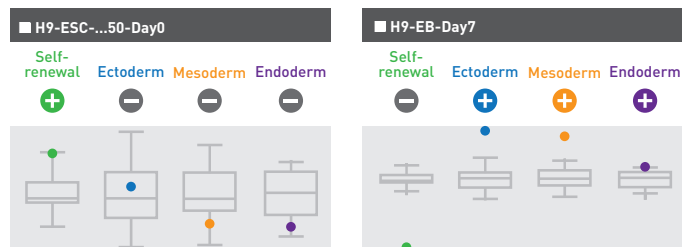
TaqMan hPSC Scorecard Panel and Analysis Software

Product	Size	Cat. No.
TaqMan™ hPSC Scorecard™ Kit, Fast 96-well	2 x 96 well plates	A15871
TaqMan™ hPSC Scorecard™ Panel, Fast 96-well	2 x 96 well plates	A15876
TaqMan™ hPSC Scorecard™ Kit, 384-well	1 x 384 well plates	A15872
TaqMan™ hPSC Scorecard™ Panel, 384-well	1 x 384 well plates	A15870

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

Product	Size	Cat. No.
PSC Neural Induction Medium	500 ml	A1647801
PSC Cardiomyocyte Differentiation Kit	1 kit	A2921201
PSC Definitive Endoderm Induction Kit	1 kit	A3062601
PSC Dopaminergic Neuron Differentiation Kit	1 kit	A3147701

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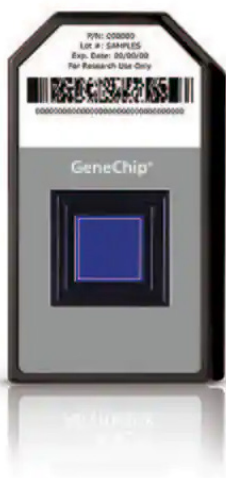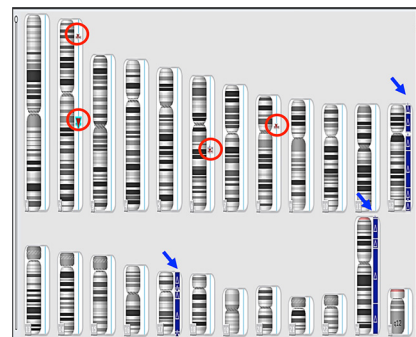Analytical Features	KaryoStat™	KaryoStat HD™
Gains	2Mb	<25-50 kb
Losses	1Mb	<25-50 kb
AOH	>5Mb	>1Mb
Mosaicism	>~20%(>30%)	>~15%(>30%)
SNP	Yes	Yes
Coverage	Low	Broad

Product	Quantity	Cat. No.
KaryoStat™ Assay	24 assays	905403
KaryoStat™ HD Assay	24 assays	905404
KaryoStat™ HD Assay Plus	24 assays	905405

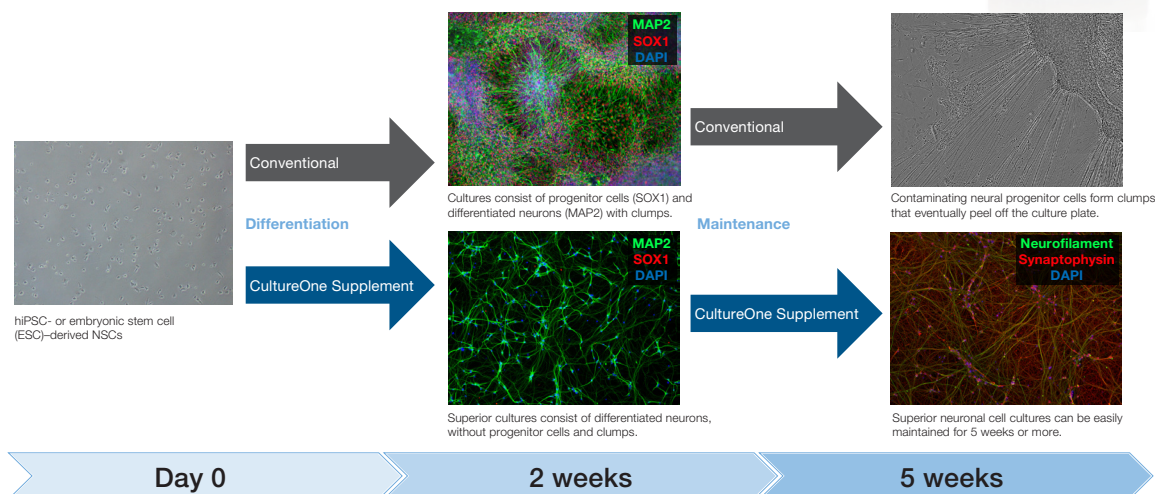
Find out more at thermofisher.com/karyostat

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
Gibco B-27™ Plus Neuronal Culture System

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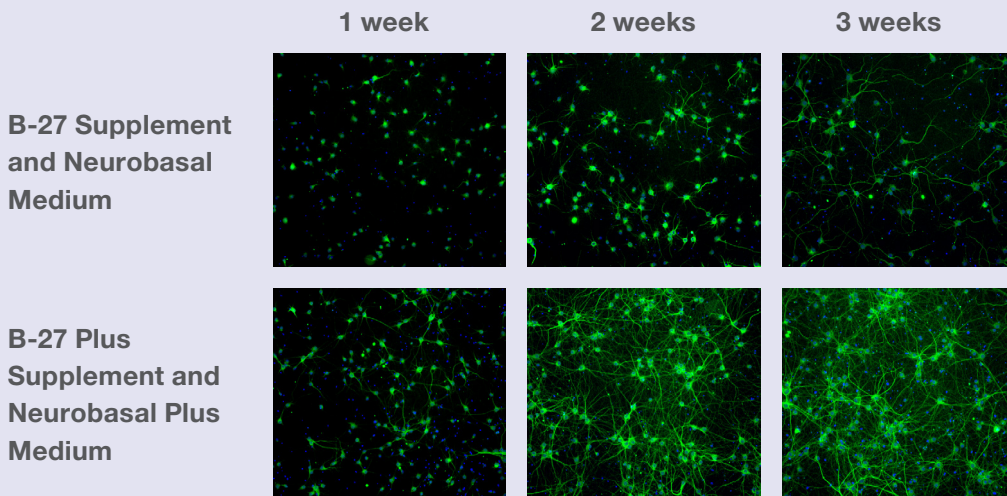- Increased neuronal survival by more than 50%
 - the highest in the industry
- Accelerated neurite outgrowth
- Improved electrophysiological activity and maturation of neurons


Product	Size	Cat. No.
B-27™ Plus Neuronal Culture System	1 system	A3653401
Neurobasal™ Plus Medium	500 ml	A3582901
B-27™ Plus Supplement (50X)	10 ml	A3582801

Learn more at thermofisher.com/b27plus

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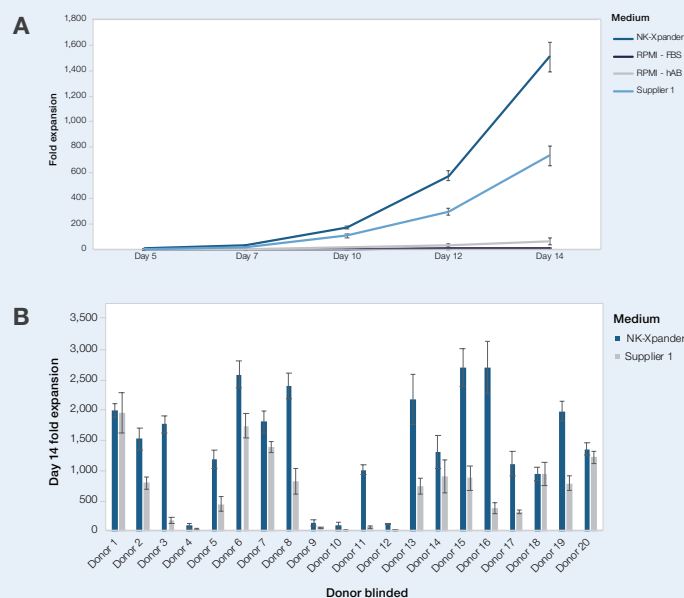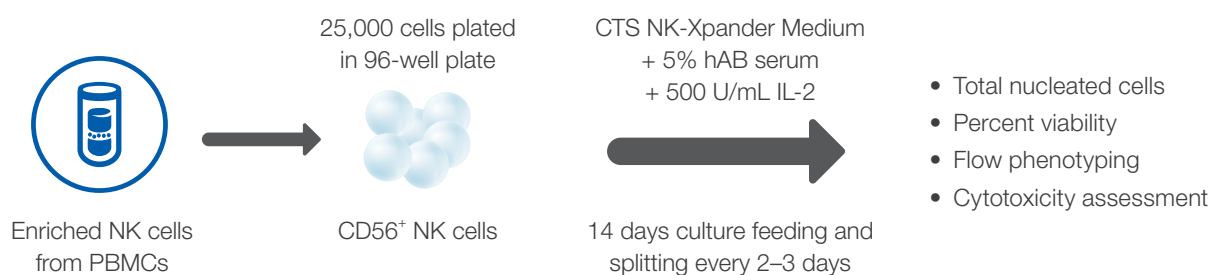
Product	Size	Cat. No.
CTST™ DPBS, calcium, magnesium	1,000 ml	A1285801

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	Cost-effective siRNA	Good knock-down, low off-target	Highest knock-down, lowest off-target
	Silencer® siRNA	Stealth RNAi siRNA	Silencer® Select siRNA
Potency	100nM recommended conc.	20nM recommended conc.	5nM recommended conc.
Guarantee	2 of 3 siRNA guaranteed	2 of 3 siRNA guaranteed	2 of 3 siRNA guaranteed
Target specificity	Moderate	High	Highest
Innate immune response	Low	Minimized through chemical modifications	Minimized through chemical modifications
Molecular format	Unmodified 21-bp duplex with overhangs	Modified 25-bp duplex with no overhangs	LNA modified 21-bp duplex with overhangs
Coverage	Coding RNA	Coding RNA	Coding & non-coding RNA
Target species	Human, mouse, rat (Other species: use custom tool)	Human, mouse, rat (Other species: use custom tool)	Human, mouse, rat (Other species: use custom tool)
Custom tool	Custom Silencer siRNA	Custom Stealth RNAi siRNA	Custom Silencer Select siRNA
Recommended Cat. No.	AM16708: Silencer Pre-designed siRNA 5nmol	1299003: RNA-Stealth Oligos Tube Set	4392420: Silencer Select Pre-Designed siRNA 5nmol

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** Silencer® Select siRNAs for noncoding RNA (ncRNA) now enable ncRNA researchers for the first time to easily obtain siRNAs targeting any ncRNA. Custom siRNA libraries for noncoding RNAs (including for lncRNAs) are available. For information on all our predefined and custom siRNA libraries, contact maillibraries@thermofisher.com.

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Learn more at thermofisher.com/sirna

Cell engineering applications and workflow solutions

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Learn more at thermofisher.com/truerecut

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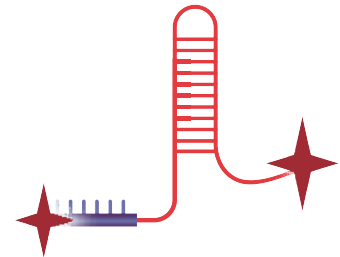
Learn more at thermofisher.com/truereguide

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Learn more at thermofisher.com/truetag

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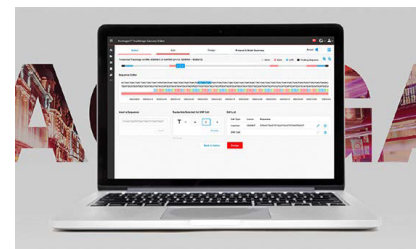
Learn more at thermofisher.com/truedesign

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
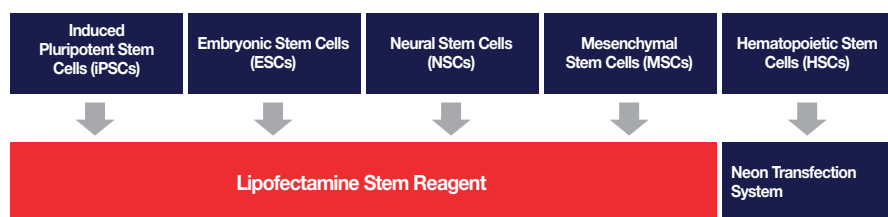
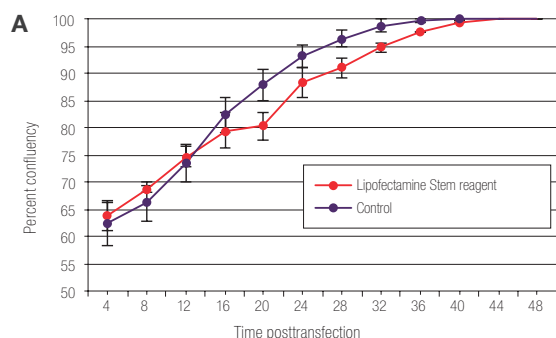
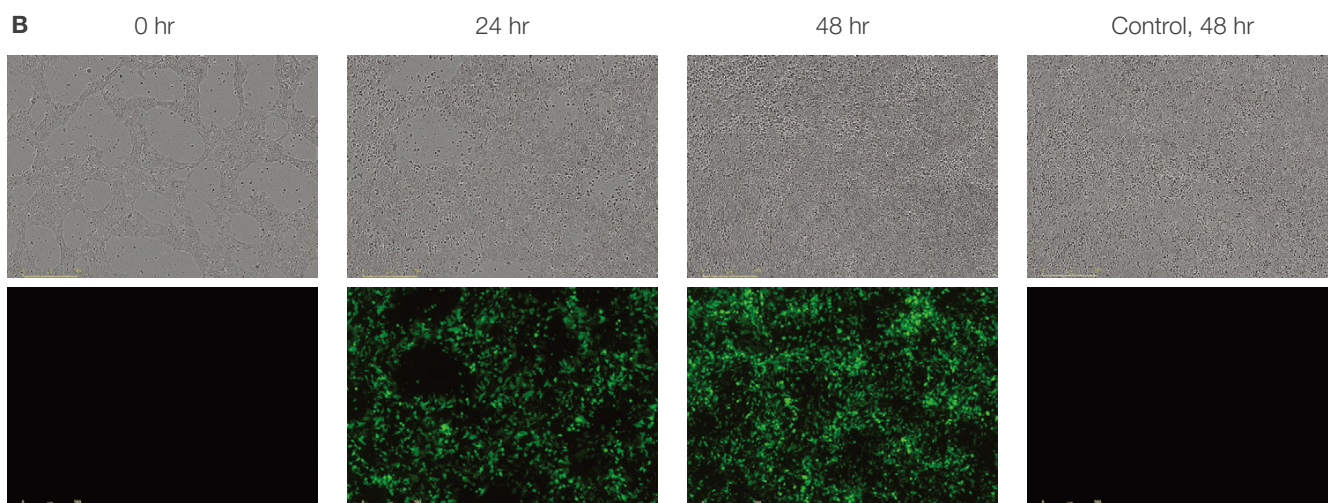
A iPSCs


Experimental condition	Recommendation
Delivery platform	Lipofectamine Stem reagent, 1 μ L/well
Plate format	24-well plate
DNA	GFP plasmid, 500 ng/well
Medium	Essential 8 Medium
Extracellular matrix	Vitronectin
Cell density	50,000 cells/well

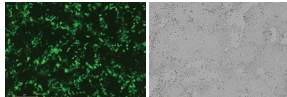iPSCs, GFP plasmid
Transfection efficiency: 75%

C iPSC-derived NSCs


Experimental condition	Recommendation
Delivery platform	Lipofectamine Stem reagent, 1 μ L/well
Plate format	24-well plate
DNA	GFP plasmid, 500 ng/well
Medium	StemPro NSC SFM
Extracellular matrix	Geltrex matrix
Cell density	75,000 cells/well

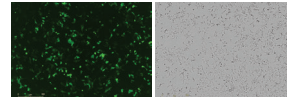NSCs, GFP plasmid
Transfection efficiency: 60%

B ESCs


Experimental condition	Recommendation
Delivery platform	Lipofectamine Stem reagent, 2 μ L/well
Plate format	24-well plate
DNA	GFP plasmid, 500 ng/well
Medium	Essential 8 Medium
Extracellular matrix	Vitronectin
Cell density	100,000 cells/well

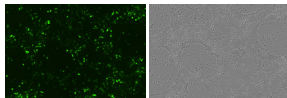H9 ESCs, GFP plasmid
Transfection efficiency: 83%


D MSCs

Experimental condition	Recommendation
Delivery platform	Lipofectamine Stem reagent, 1 μ L/well
Plate format	24-well plate
DNA	GFP plasmid, 500 ng/well
Medium	MesenPRO RS Medium
Extracellular matrix	CTS CELLstart Substrate
Cell density	25,000 cells/well

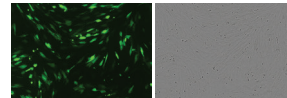Adipose-derived MSCs,
GFP plasmid
Transfection efficiency: 47%

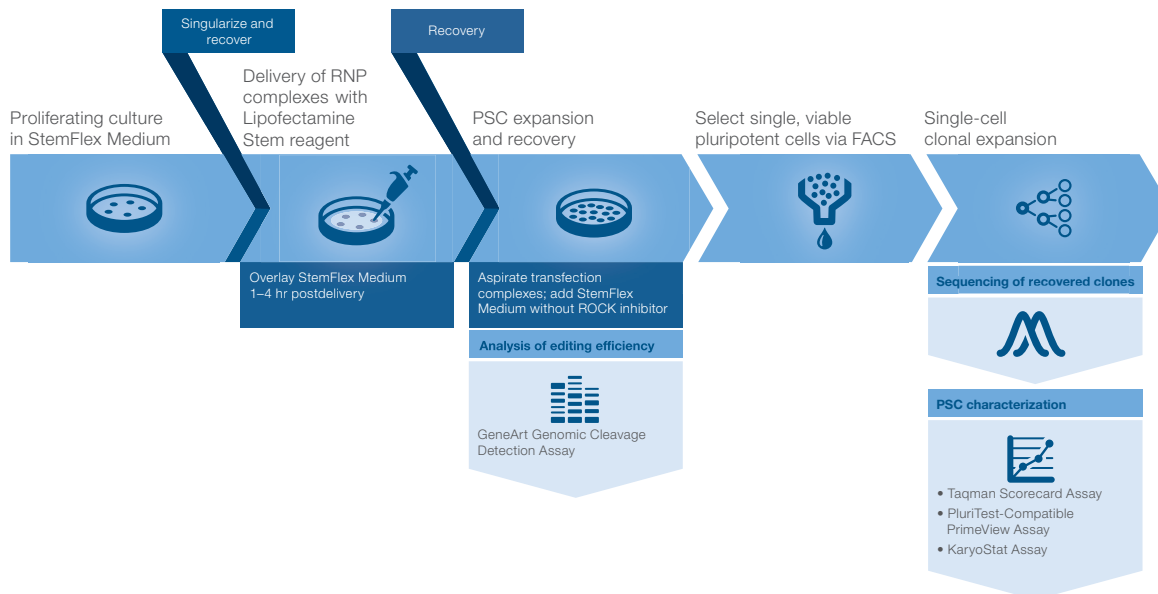Invitrogen™ Lipofectamine™ Stem Transfection Reagent의 gene editing application workflow

Product	Size	Cat. No.
Lipofectamine™ Stem Transfection Reagent	0.1 ml	STEM00001
	0.3 ml	STEM00003
	0.75 ml	STEM00008
	1.5 ml	STEM00015

CTS™ Xenon Electroporation System

Closed, optimized, and scalable platform for electroporation

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
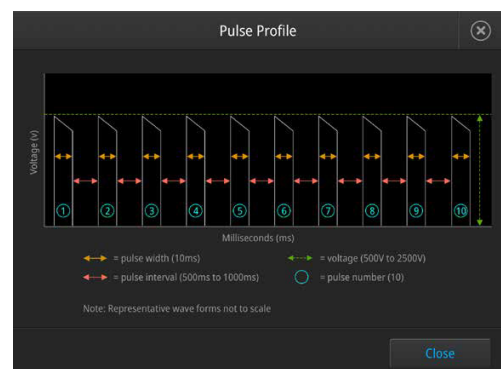
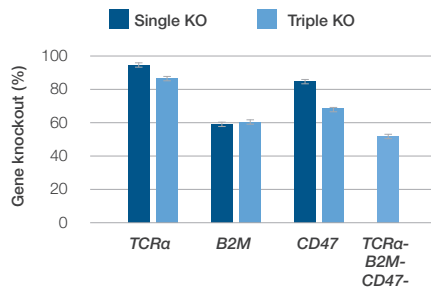
Large scale Xenon Electroporation System

Seamless translation

- Day 3 T cells (5×10^7 cells/mL)
- Analysis performed 72 hr post-EP
- Single KO payload mix :
- Cas9 v2 : 120 µg/mL
- gRNA : 28 µg/mL
- Triple KO payload mix :
- Cas9 v2 : 360 µg/mL
- gRNA : 28 µg per target
- Genome Editing Buffer

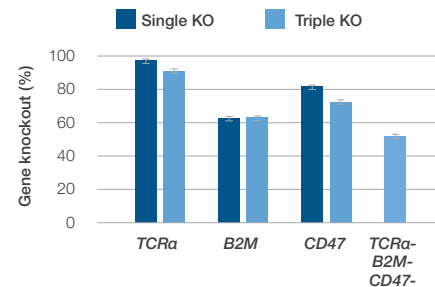Small scale Neon Transfection System

Recommended process volume cell concentration	20×10^6 to 100×10^6
Electroporation volumes	1 mL (CTS™ Xenon SingleShot chamber) 5–25 mL (Xenon MultiShot cartridge)
Electroporation chamber volume	1 mL
Run time for 5–25 mL electroporation volume	7–22 min
Electroporation pulse voltage range	500–2,500 V
Electroporation pulse width range	1–30 ms
Electroporation pulse interval range	500–1,000 ms
Number of electroporation pulses	1–10
Cell mixer rotation speed	60 rpm
Precooling technology	Peltier
Precooling temperature setting range	10–30°C

 Learn more at [thermofisher.com](https://www.thermofisher.com)

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For Research Use Only. Not for use in diagnostic procedures. © 2022 Thermo Fisher Scientific Inc.
All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

ThermoFisher
S C I E N T I F I C