WESTERN DETECTION

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The Invitrogen™ iBlot™ 3 Western Blot Transfer System is an essential part of the Invitrogen™ western blotting workflow.

Electrophoresis

Protein gels welcome packs

Invitrogen™ PowerEase™ Touch power supplies

Transfer

iBlot 3 Western Blot Transfer System

Processing

Invitrogen™ iBind™ and iBind™ Flex automated western processing devices

Primary and secondary antibodies

Imaging

Invitrogen™ iBright™ Imaging Systems

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Figure 1. The all-new design of the iBlot 3 device features two independently controlled transfer stations with programmable cooling.

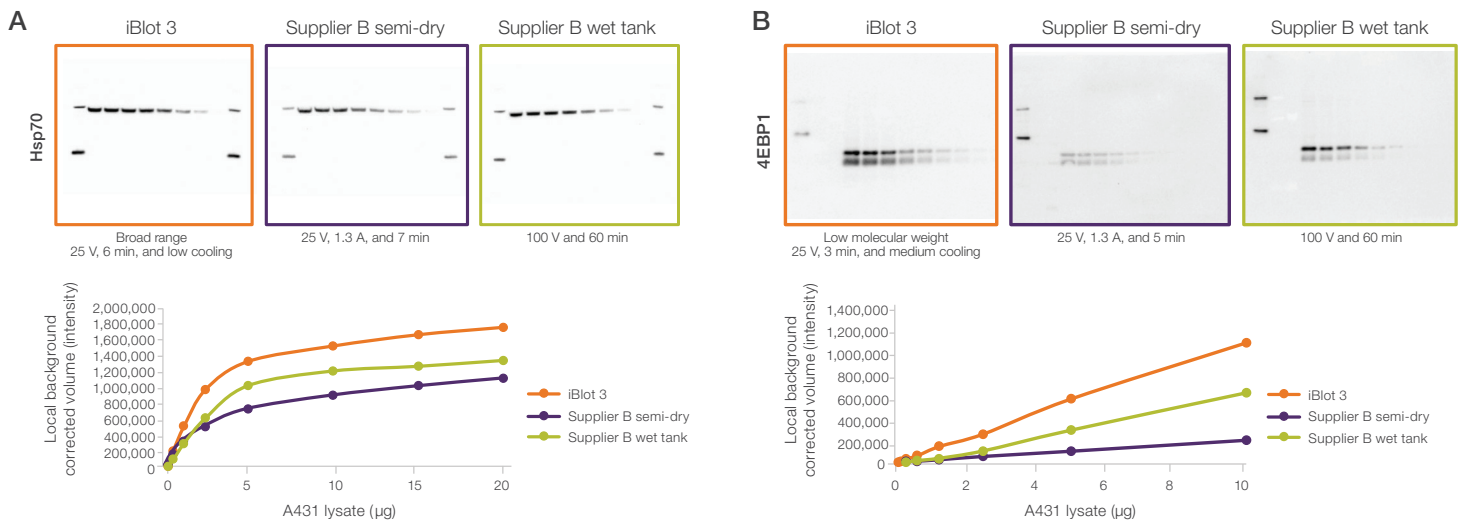
Figure 2. Four mini transfer stacks loaded in the iBlot 3 device.

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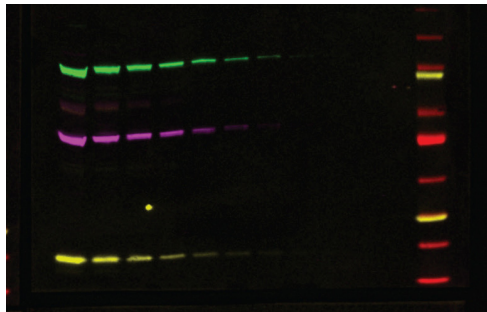
Figure 4. Invitrogen™ iBolt™ 3 Transfer Stacks.
Each box contains 10 stacks (mini or midi format),
10 filter papers, and 10 absorbent pads.

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Transfer station 1 with two gels

Transfer station 2 with two gels

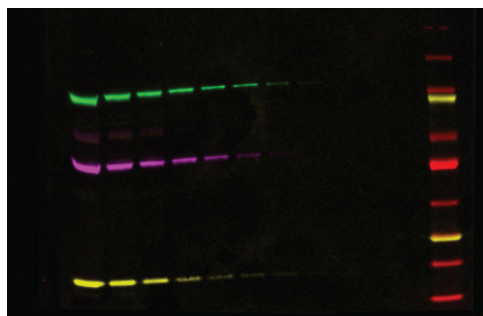
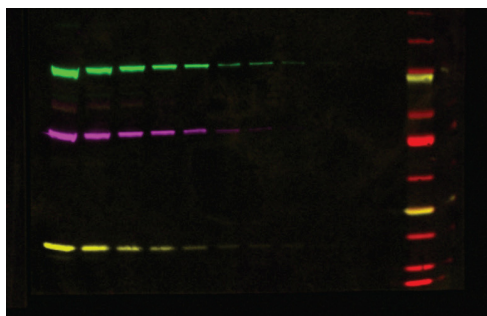
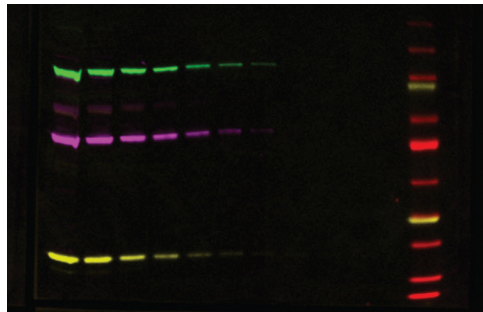
Figure 5. The iBolt 3 device delivers consistent results. A serial dilution of HEK293 lysate (lysed in Thermo Scientific™ RIPA buffer) was loaded onto four Invitrogen™ Bolt™ 4 to 12% Bis-Tris Plus gels. Protein loads were (left to right, in μg): 10, 5, 2.5, 1.25, 0.625, 0.312, 0.156, 0.078, 0.039, 0.020, and 0.010. Invitrogen™ iBright™ Prestained Protein Ladder was loaded into the last lane. MES-SDS running buffer was used for gel electrophoresis. Proteins were transferred using the iBolt 3 Western Blot Transfer Device and iBolt 3 Transfer Stacks, Midi, NC using the broad range (30–250 kDa) preprogrammed transfer method (25 V, 6 min, low cooling). Blots were then blocked with Pierce Clear Milk Blocking Buffer and were then probed using primary antibodies raised against epidermal growth factor receptor (EGFR), calreticulin, and p23 in clear milk for 12 hours. Secondary antibodies Invitrogen™ GAR-Alexa Fluor™ Plus 800 dye, Invitrogen™ GAM-Alexa Fluor™ Plus 488 dye, and Invitrogen™ GACHk-Alexa Fluor™ 546 dye were prepared in TBST and added to the blots for 2 hours. Immunoprocessing was completed on the Bandmate Automated Western Blot Processor. Imaging was completed on the iBright FL1500 Imaging System. Note: GACHk = goat anti-chicken.

Ordering information

Description	Quantity	Cat. No.
iBlot 3 Western Blot Transfer Device	1	IB31001
iBlot 3 Transfer Stacks, Midi, NC	10 stacks	IB33001
iBlot 3 Transfer Stacks, Mini, NC	10 stacks	IB33002
iBlot 3 Transfer Stacks, Midi, PVDF	10 stacks	IB34001
iBlot 3 Transfer Stacks, Mini, PVDF	10 stacks	IB34002
iBlot 3 Starter Kit, NC	1 kit	IB31001S
iBlot 3 Starter Kit, PVDF	1 kit	IB31002S
iBlot 3 IBind/iBind Flex Starter Kit*	1 kit	IB31001SLF1/ IB31001SLF2
Bolt Welcome Pack with iBlot 3 system	1 pack	NW0412AIB3
Starter kit and welcome pack contents		
iBlot 3 Starter Kit, NC		
iBlot 3 Western Blot Transfer Device	1	IB31001
iBlot 3 Transfer Stacks, Midi, NC	2 boxes (20 transfer stacks)	IB33001
iBlot 3 Starter Kit, PVDF		
iBlot 3 Western Blot Transfer Device	1	IB31001
iBlot 3 Transfer Stacks, Midi, PVDF	2 boxes (20 transfer stacks)	IB34001
iBlot 3 IBind Starter Kit		
iBlot 3 Western Blot Transfer Device	1	IB31001
iBlot 3 Transfer Stacks, Midi, NC	1 box (10 transfer stacks)	IB33001
iBind Western Device	1	SLF1000
iBind Cards	1	SLF1010
iBind Solution Kit	1	SLF1020
SuperSignal West Pico Plus Chemiluminescent Substrate	200 mL	34577
Bolt Welcome Pack with iBlot 3 system		
iBlot 3 Western Blot Transfer Device	1	IB31001
iBlot 3 Transfer Stacks, Midi, NC	1 box (10 transfer stacks)	IB33001
Mini Gel Tank	1	A25977
20X Bolt MES SDS Running Buffer	500 mL	B0002
4X Bolt LDS Sample Buffer	10 mL	B0007
10X Bolt Sample Reducing Agent	10 mL	B0009
PageRuler Plus Prestained Protein Ladder, 10–250 kDa	2 x 250 µL	26619
Bolt 4 to 12% Bis-Tris 1.0 mm, Mini Protein Gels	1 box (10 gels)	NW04120BOX

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