Simply brilliant

One-touch gel imaging

iGlow Gel Documentation System for Protein and Nucleic Acid Gels

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invitrogen

iGlow Gel Documentation System— one-touch gel imaging that's simply brilliant

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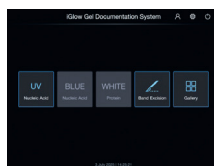
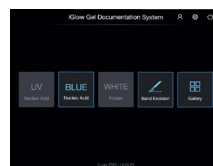
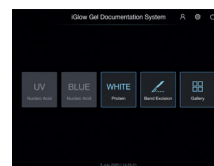
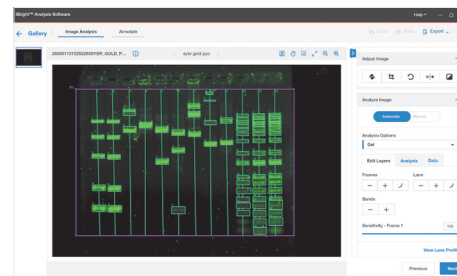
iGlow system stains and dyes compatibility chart

Application	Gel stain	Excitation source		
		UV	Blue*	White
Protein	Invitrogen™ Coomassie Fluor™ Orange		X	
	BioRad™ Flamingo™ Stain		X	
	BioRad™ Oriole™ Stain	X		
	Thermo Scientific™ Pierce™ Silver Stain			X
	Invitrogen™ Pro-Q™ Diamond Phosphoprotein		X	
	Invitrogen™ SimplyBlue™ SafeStain			X
	Invitrogen™ SYPRO™ Ruby	X		
	Invitrogen™ SYPRO™ Orange	X		
	Invitrogen™ SYPRO™ Red	X		
	Invitrogen™ SYPRO™ Tangerine	X		
DNA	Millipore-Sigma™ GelRed™ stain	X		
	Gold-Bio™ GelGreen™ stain		X	
	Bulldog Bio™ Midori Green		X	
	Invitrogen™ SYBR™ Gold		X	
	Invitrogen™ SYBR™ Green		X	
	Invitrogen™ SYBR™ Safe		X	
	Invitrogen™ UltraPure™ Ethidium Bromide	X		

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
Learn more at thermofisher.com/iglow

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
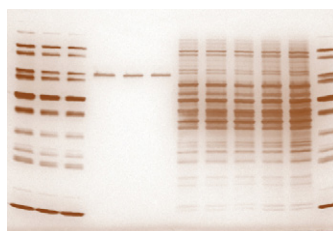
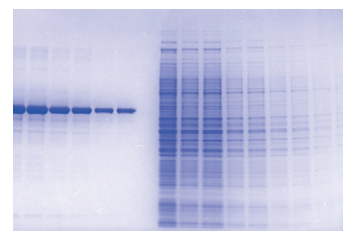
Stain used: Pierce Silver Stain (Cat. No. 24612)
Light source: white

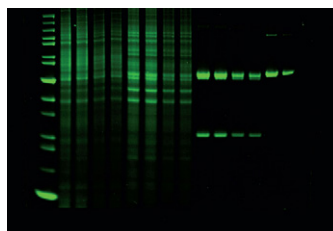Stain used: SimplyBlue SafeStain (Cat. No. LC6065)
Light source: white

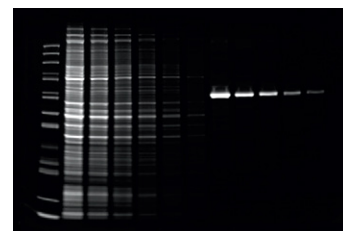
Stain used: SYPRO Red Protein Gel Stain (Cat. No. S6653)
Light source: UV


Stain used: Coomassie Fluor Orange (Cat. No. C33250)
Light source: blue

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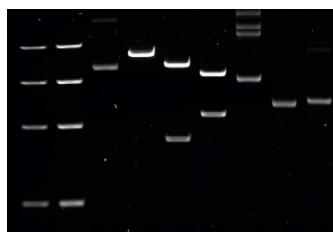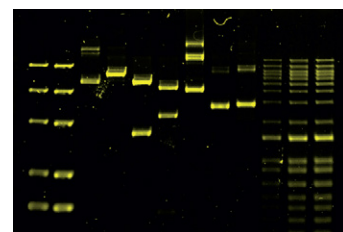Stain used: UltraPure Ethidium Bromide (Cat. No. 15585011)
Light source: UV

Stain used: SYBR Gold Nucleic Acid Gel Stain (Cat. No. S11494)
Light source: blue

Technical specifications

Optics	
Imaging modes	Colorimetric, fluorescent, GFP colony counting
Detector	Scientific CMOS (sCMOS)
Resolution	5.0 megapixels
Field of view	5.5 x 8.3 in. (14 x 21 cm)
System interface	
Touchscreen	12.1 in. LCD display, 1,024 x 768 pixels
Sample drawer	Manually operated
Connectivity and storage	
USB version	2.0, 3.0
Networking	Ethernet port, 2.4 GHz Wi-Fi
Image file types	TIFF, JPG, PNG, IBA (proprietary)
Built-in computer details	
Hard drive	500 GB solid-state drive
Storage available	400 GB
RAM	8 GB
Illumination sources	
Transillumination	UV, blue, and white LED
LED spectra	UV: 300–310 nm, Blue: 460–470 nm, White: 400–800 nm
Preinstalled filter	590 nm emission filter
Physical characteristics	
Power supply	100–240 VAC
Dimensions (L x W x H)	15 x 18 x 16 in. (38 x 45 x 41 cm)
Weight	48.5 lb (22 kg)
Operating conditions (indoor use)	5–40°C, <80% humidity
Support	
Standard warranty	2 years from date of purchase

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For Research Use Only. Not for use in diagnostic procedures. © 2026 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. **SPEC-13605369 0326**