PCR

New products across the PCR workflow

Invitrogen™ SuperScript™ Reverse Transcriptases

Invitrogen™ Platinum™ DNA Polymerases

Applied Biosystems™ Thermal Cycler

15%

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STEP 1

cDNA Synthesis

Which SuperScript IV format is right for you?

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SuperScript IV First-Strand Synthesis System

- 합성에 필요한 모든 **Component**가 포함
- 추가적으로 Hela RNA, Control gene primer set 제공
- 구성품 : SuperScript® IV RT, 5X RT buffer, 0.1 M DTT, 10 mM dNTP mix, Oligo(dT)20, Random hexamers, Ribonuclease Inhibitor, E. coli RNase H, DEPC-treated water, Total HeLa RNA, Control Primer set

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SuperScript IV One-Step RT-PCR System

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Why should you choose SuperScript IV Reverse Transcriptases?

SuperScript IV Reverse Transcriptase (RT) features:

Super Fast

Short, 10 min cDNA synthesis protocol

Super Efficient

Higher cDNA yields than with other RT enzymes

Super Stable

High thermostability and processivity for superior cDNA synthesis performance

Super Strong

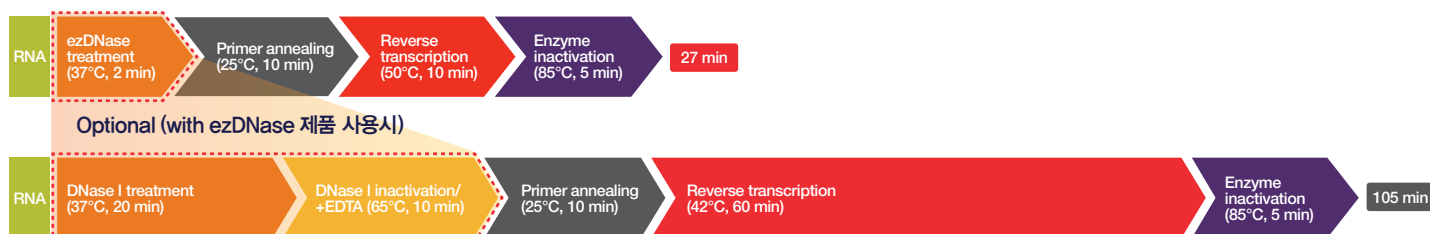
Great results, even with RNA samples of suboptimal purity

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Traditional cDNA synthesis workflow with DNase I

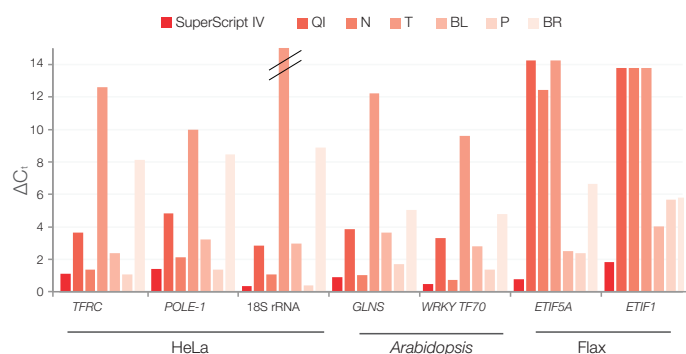
Figure 1. High efficiency with degraded RNA. RT-qPCR of degraded RNA (RNA integrity number (RIN) 1–3) from human cells and plant tissues with different RTs and Applied Biosystems™ TaqMan® Assays. Delta C_t values ($\Delta C_t = C_t - C_{t \text{ SuperScript IV}}$) show that SuperScript IV RT, with its standard 10 min protocol, delivered higher cDNA yields and lower C_t values than the recommended protocols for SuperScript III reagent and other suppliers' RTs.

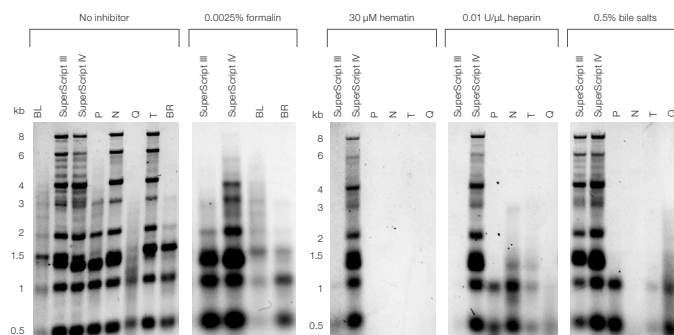
Figure 2. Tolerance to inhibitors. This data shows first-strand cDNA synthesis of an RNA ladder in the presence of reaction inhibitors. cDNA fragments were resolved by alkaline gel electrophoresis. All RTs except for SuperScript IV RT were severely affected with trace amounts of inhibitors.

STEP 2

DNA Amplification

Performance and flexibility to fit your amplification

Unique features of the latest Platinum DNA polymerases

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

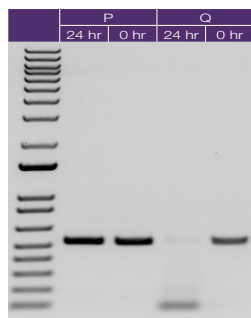
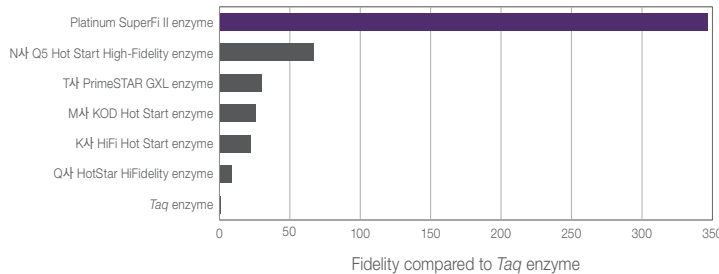
Figure 3. Fidelity comparison of commercially available enzymes and Taq enzyme. PCR amplicons (3.9 kb) obtained using different DNA polymerases were fragmented with a MuA transposase. Unique molecular identifiers (UMIs) of 12 random nucleotides were introduced during fragmentation to tag each product individually. After next-generation sequencing, reads from the same UMI family were aligned to call errors. Errors were identified only when present in all reads in the UMI family; otherwise they were discarded as sequencing errors. The fidelity values were normalized to Taq polymerase fidelity.

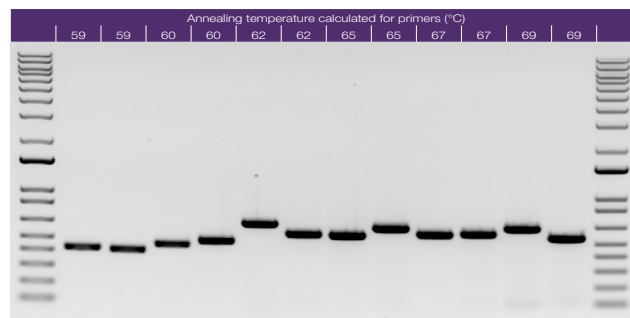Benchtop stability for high-throughput applications

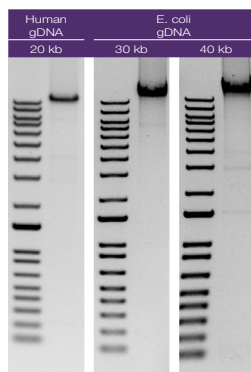
No need to calculate primer annealing temperature

High success with amplification of 20–40 kb sequences

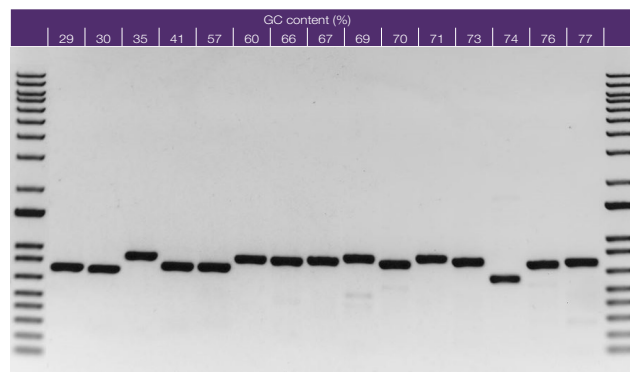Figure 4. Amplification of long fragments. A 20 kb target from human gDNA and 30–40 kb targets from E. coli gDNA were successfully amplified using Platinum SuperFi II DNA Polymerase.

No additives needed for GC-rich amplification

Figure 5. Robust amplification of GC-rich targets. Fifteen targets with a range of GC content were amplified from human gDNA without any supplementary buffer additives.

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Fast DNA synthesis at 15 sec/kb

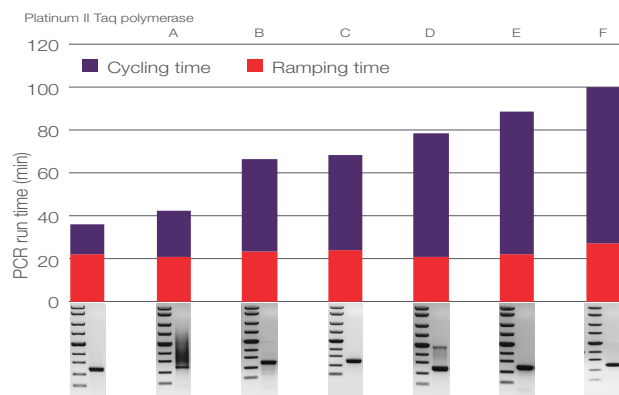
Figure 6. Fast cycling reduces PCR run time. A 0.5 kb fragment was amplified for 35 cycles, using Platinum II Taq Hot-Start DNA Polymerase and hot-start DNA polymerases from other suppliers (A–F).

High sensitivity and specificity

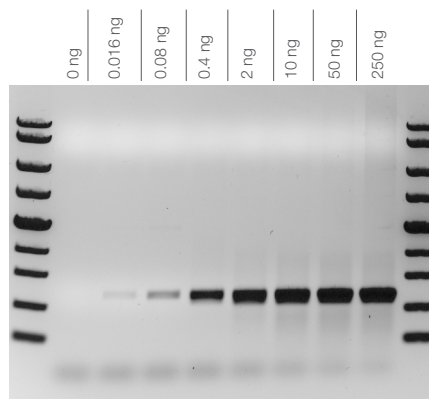
Figure 7. High sensitivity and reliable amplification from low amounts of input DNA. A 0.5 kb fragment was amplified from a range of human gDNA input, using Platinum II Taq Hot-Start DNA Polymerase.

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Two protocols to fit your needs

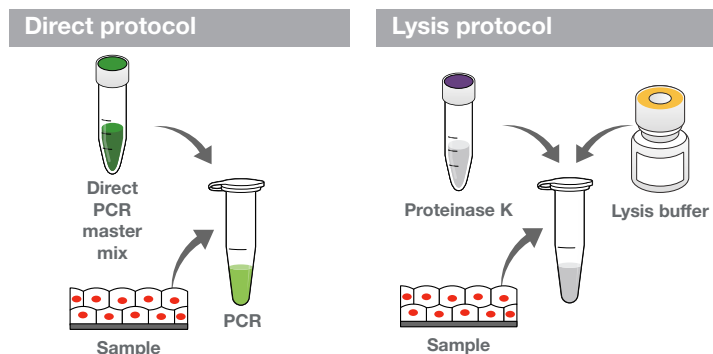
Figure 8. Direct and lysis protocols. Two protocols are available to amplify target DNA directly from crude samples. The direct protocol offers a shorter workflow, whereas the lysis protocol allows flexibility and long-term storage.

Multiple PCR targets in the same reaction

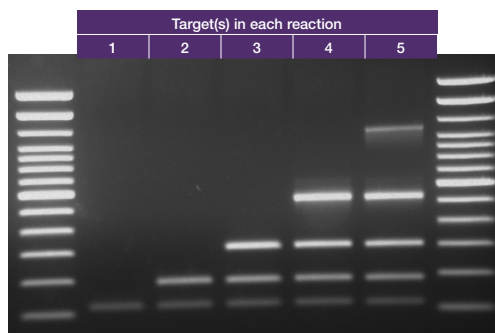
Figure 9. Efficient multiplexing in direct PCR. Five fragments of 0.1–1.1 kb were amplified from mouse tail in singleplex to 5-plex reactions, using the lysis protocol.

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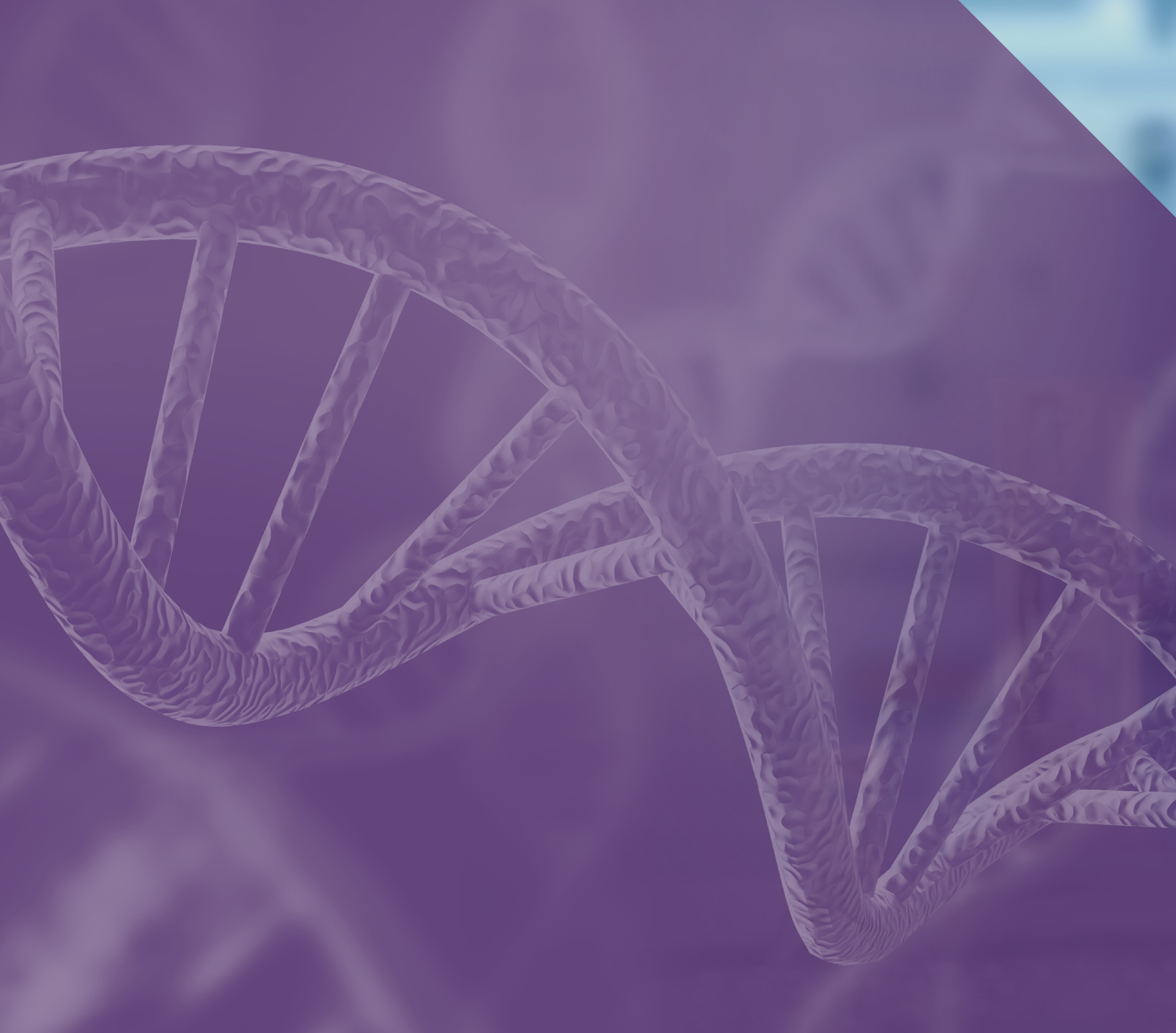
For Research Use Only. Not for use in diagnostic procedures. © 2021 Thermo Fisher Scientific Inc.
All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

ThermoFisher
S C I E N T I F I C